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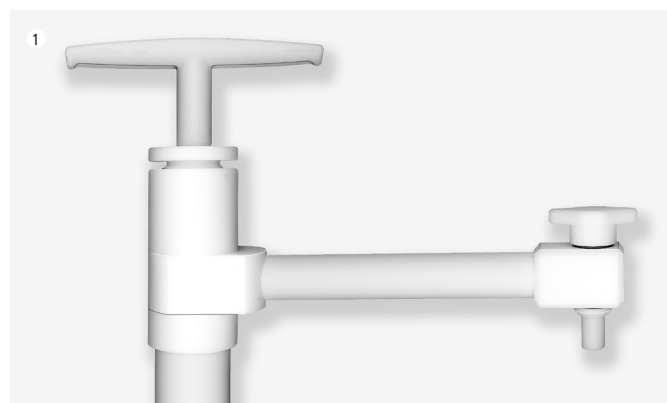
# Instructions for Use PTFE Barrel pump ultraclean

Art. No. 5606-1000 | 5606-0600 Read carefully before use and keep for later reference.

## Safety regulations

- ▶ When transferring flammable liquids please observe the "Technical Rules for Flammable Liquids" (TRBF) and the safety hints given by the reagent manufacturer (safety data sheet).
- ▶ Protective clothing must always be worn handling hazardous liquids (protective goggles, gloves, apron and, if necessary, boots).
- ▶ Each user must know about these instructions and have them to hand at all times.
- ▶ Only use the pump in accordance with its resistance properties. Medium-contacting parts are made from PTFE and FEP.
- ▶ Gas-emitting and fuming media should only be transferred under fume hoods or in well-ventilated rooms. Observe the safety hints given by the reagent manufacturer (safety data sheet).
- ▶ Always handle the apparatus so that neither the user nor other persons are at risk.
- ▶ Never use force.

## Packaging/Contents

Check the contents for completeness and any transport damage. If external damage to the packaging is obvious then immediately inform the deliverer/ forwarding agent and obtain confirmation.

**Warning: Avoid injury by hazardous liquids! Parts of the apparatus could still contain the liquid. Never point openings toward your body! Wear protective goggles, gloves and apron! Do not overfill the receiving container!**

## Function, safety measures, limitations of use

This pump is used to transfer low-viscosity to oily liquids from canisters and drums up to approx. 220 liters. It is the ideal solution for transferring high-purity liquids. The universal chemical resistance of the medium-contacting materials PTFE and FEP means that liquids which attack conventional materials such as V2A steel, PP or PVC can be transferred.

The stopcock stops the liquid from dripping as well as the escape of hazardous vapors. This why the stopcock should only be opened during the transfer process, see Fig. 3. When the transfer process is finished the stopcock should be returned to the closed position, see Fig. 2.

Keep the container to be filled away from your body and wear protective clothing. The user must check that the apparatus is suitable for the particular application.

After you have finished filling, push piston rod into the pump completely.

## Cleaning and maintenance

Please take care that the liquid does not contain any large particles of dirt, as these could destroy the pump valves and passages. Rinse the pump with a suitable medium at regular intervals! If the pump is damaged it should no longer be used and returned for repair!

## Accessory

	Art-No.
Full-protection goggles, non misting	2502-3001
Protective gloves, nitrile	2507-0004
Protective apron	2507-1001
Safety stand	5618-0100
Universal barrel wrench	5690-0090
Antistatics-Set	5602-1000
Barrel screw joint nickel-plated brass R2" internal fine thread, steel	5601-0100
Barrel screw thread, gas tight	5601-0130
Barrel screw joints, PP Mauser 2", internal coarse thread	5600-3170
Tri-Sure, internal coarse thread	5600-3180
PVC Barrel stopper for 40-70 mm dia. openings	5600-3150

## Faults – troubleshooting

Liquid is not transferred:

- base valve dirty

If the container contains only a little residual liquid then pumping must be carried out for a longer period.